

# ETL500 Tube holder setting guide

## ETL500 pumphead - Instruction for tube holders



Open the pump by moving the lever (flip top) to an upright position.

Set tube holders on both sides by turning the adjustment wheels.

See tables below for recommended positions for standard tube material and inner diameter sizes.

### Pharm-A-Line tube (TPV)

Tube bore (ID mm)	Scale position
0.5	1
0.8	1
1.6	2
3.2	3
4.8	4
6.4	6
8.0	7



### Pharm-A-Clear tube (Silicone)

Tube bore (ID mm)	Scale position
0.5	1
0.8	1
1.6	2
3.2	3
4.8	4
6.4	6
8.0	7



Concerning other tube materials and/or sizes then the above the scale position the tube holder setting needs to be tested out, but present scale positions can be used as first input for a functional test. Wall thickness of the tube needs to be 1.6mm for this pump version

**THOMAS**  
An Ingersoll Rand Business

[www.thomaspumps.com](http://www.thomaspumps.com)